# RotoForce-3

Instruction Manual

Manual No.: 15247000

Date of Release 30.01.2001

*RotoForce-3*  
*Instruction Manual*

<b>Table of Contents</b>	<b>Page</b>
<b>User's Guide</b> .....	<b>1</b>
<b>Reference Guide</b> .....	<b>15</b>
<b>Quick Reference Guide</b> .....	<b>23</b>

---

Always state *Serial No* and *Voltage/frequency* if you have technical questions or when ordering spare parts. You will find the *Serial No.* and *Voltage* on the type plate of the machine itself. We may also need the *Date* and *Article No* of the manual. This information is found on the front cover.

The following restrictions should be observed, as violation of the restrictions may cause cancellation of Struers legal obligations:

**Instruction Manuals:** Struers Instruction Manual may only be used in connection with Struers equipment covered by the Instruction Manual.

**Service Manuals:** Struers Service Manual may only be used by a trained technician authorised by Struers. The Service Manual may only be used in connection with Struers equipment covered by the Service Manual.

Struers assumes no responsibility for errors in the manual text/illustrations. The information in this manual is subject to changes without notice. The manual may mention accessories or parts not included in the present version of the equipment.

The contents of this manual is the property of Struers. Reproduction of any part of this manual without the written permission of Struers is not allowed.

All rights reserved. © Struers 2001.

**Struers A/S**  
Valhøjs Allé 176  
DK-2610 Rødovre/Copenhagen  
Denmark  
Telephone +45 36 70 35 00  
Fax +45 38 27 27 01

---

## **RotoForce-3**

### **Safety Precaution Sheet**

#### **To be read carefully before use**

1. The operator should be fully aware of the use of the machine according to the Instruction Manual. The machine must be placed in an adequate working position.
2. Be sure that the actual voltage corresponds to the voltage stated on the back of the machine. The machine must be earthed.
3. The specimen holder disc must be clamped in the quick coupling prior to starting the machine. RotoForce-3 must be correctly and securely mounted on the grinding/polishing machine.
4. Do not touch the rotating parts during operation.
5. If you observe malfunctions or hear unusual noises - stop the machine and call technical service.
6. Blue lubricant: follow the current safety rules for handling, mixing, filling, emptying and disposal of the alcohol-based lubricant.

---

The equipment is designed for use with consumables supplied by Struers. If subjected to misuse, improper installation, alteration, neglect, accident or improper repair, Struers will accept no responsibility for damage(s) to the user or the equipment.

Dismantling of any part of the equipment, during service or repair, should always be performed by a qualified technician (electromechanical, electronic, mechanical, pneumatic, etc.).

---

# User's Guide

Table of Contents	Page
<b>1. Getting Started</b>	
Checking the Contents of Packing .....	2
Unpacking of RotoForce-3 .....	2
Getting Acquainted with RotoForce-3.....	2
Installation.....	3
RotoForce-3 on RotoPol-31.....	3
RotoForce-3 on RotoPol-21/22/25 or on Rotopol-1/-2/-V .....	3
Connecting RotoForce-3.....	4
Electrical Connection.....	4
Connection to Compressed Air .....	4
Connection to the Preparation Equipment .....	4
Choosing between Menus .....	5
Setting the Language.....	6
Mounting Lupo.....	7
<b>2. Operation</b>	
Using the Controls.....	8
Front Panel Controls of RotoForce-3 and Lupo .....	8
Front Panel Controls .....	9
RotoForce-3 .....	9
Lupo.....	9
Location of Main Switch .....	10
General Use of Keys .....	10
Acoustic Signals.....	10
The Handle.....	10
Changing to Another Lupo Unit (Accessory) .....	10
Changing Lubricant Type / Refilling.....	11
Running a Preparation Method.....	12
Inserting a Specimen Holder .....	12
Adjusting the Position of RotoForce-3.....	12
Adjusting the Force.....	12
Adjusting the Preparation Time.....	13
Operation of Lupo .....	13
Operation of RotoForce-3.....	13

## 1. Getting Started

### Checking the Contents of Packing

In the packing box you should find the following parts:

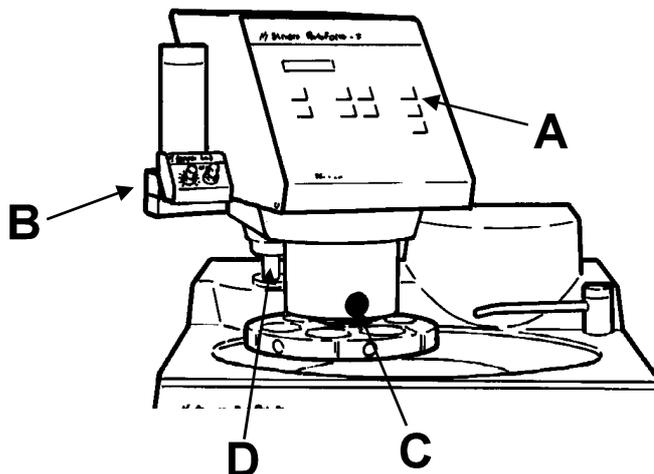
- 1 RotoForce-3
- 1 Lupo automatic drip lubricator
- 1 Silicone tube
- 1 Pressure hose, 3 m
- 2 Hose clamps
- 1 Hose clamp
- 1 Quick coupling
- 1 Hose nipple
- 1 Screw M5x10 with washer
- 1 Socket spanner
- 1 Allen key
- 1 Set of Instruction manuals

### Unpacking of RotoForce-3

Remove RotoForce-3 from the transportation box. Save the nut and the washer holding RotoForce-3. They are to be re-used when mounting RotoForce-3 on RotoPol-21/-22/-25 or Rotopol-1/-2/-V.

### Getting Acquainted with RotoForce-3

Take a moment to familiarise yourself with the location and names of the RotoForce-3 components.

*RotoForce-3 on RotoPol-31 (optional)*

- A Control keys and display
- B Lupo drip lubricator
- C Release knob
- D Supporting column and stop ring

## **Installation**

### *RotoForce-3 on RotoPol-31*

- Lift off the cap covering the supporting hole placed on the top of the RotoPol-31 cabinet with a knife.
- Likewise remove the two caps covering the fastening screws on the left-hand side of the RotoPol cabinet.
- Remove the black stop ring from the supporting column of RotoForce-3.
- Lead the supporting column of RotoForce-3 down into the supporting hole in RotoPol-31. Turn the column until the cross pin in the bottom of the column gets caught by the guiding slots inside RotoPol-31.
- Tighten the two fastening screws on the left-hand side of the RotoPol-31 cabinet and cover the screws again with the two caps.

### *RotoForce-3 on RotoPol-21/22/25 or on Rotopol-1/-2/-V*

- Place two working tables of equal height at a distance about half the length of RotoPol. Place RotoPol between the tables, carefully balancing the machine, until it stands firmly between the tables.

***Take care!***

A RotoPol machine equipped with a RotoForce weights over 50 kg!

- *RotoPol-21/-22/-25:* Remove the cap covering the supporting hole between the two turntables.  
*RotoPol-1/-2/-V:* The transportation case for RotoPol includes a printed template. Punch out the pre-cast hole in RotoPol with a round-headed hammer as follows: strike the circle on the template marked *RotoForce-3* with a firm blow. Remove the template when the hole opens.
- Lead the supporting column of RotoForce-3 down into the supporting hole of RotoPol.
- Turn the column until the cross pin in the bottom of the column gets caught by the guiding slots inside RotoPol.
- Fasten the nut and washer from the transportation box tightly under the column of RotoForce-3 below RotoPol.

## **Connecting RotoForce-3**

RotoForce-3 is supplied with electric cable, air hose and connection cable.

### *Electrical Connection*

Mount a plug on the cable or connect directly to mains:

Black and brown: phase

Yellow/green: earth

**IMPORTANT**

Check that the mains voltage corresponds to the voltage stated on the type plate on the back of the machine.

Check that the specimen holder rotates anti-clockwise.

### *Connection to Compressed Air*

- Mount an air hose on the enclosed quick coupling and secure with the enclosed hose clamp.
- Connect the quick coupling to the factory-mounted air hose of RotoForce-3.

**IMPORTANT**

The air pressure must be between 5 bar (73 psi) and 6.5 bar (95 psi)  
The air must not contain water, oil or particles. If necessary, a pressure regulator/filter unit must be installed between air supply and RotoForce.

### *Connection to the Preparation Equipment*

*RotoPol alone:* Connect the communication cable to the back of RotoPol (see the Instruction Manual of RotoPol).

*RotoPol with Multidoser and/or RotoCom:* Connect the communication cable to the back of Multidoser (see the Instruction Manual of Multidoser)

**Choosing between Menus**

Press the MENU  key to move from one screen to another:

LANGUAGE  
ENGLISH

COMMUNICATION  
STANDARD

FORCE REDUCTION  
NO

## Setting the Language

150N 01:00  
STOP

LANGUAGE  
ENGLISH

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150N 01:00  
STOP

Press MENU  to activate the set-up menu.

The LANGUAGE screen appears.

Press SELECT  to choose English, German, French or Japanese.

Press STOP  to accept the choice. The Operation Menu will appear again in the language you have chosen.

Remember to set the same language on all the connected RotoSystem modules.

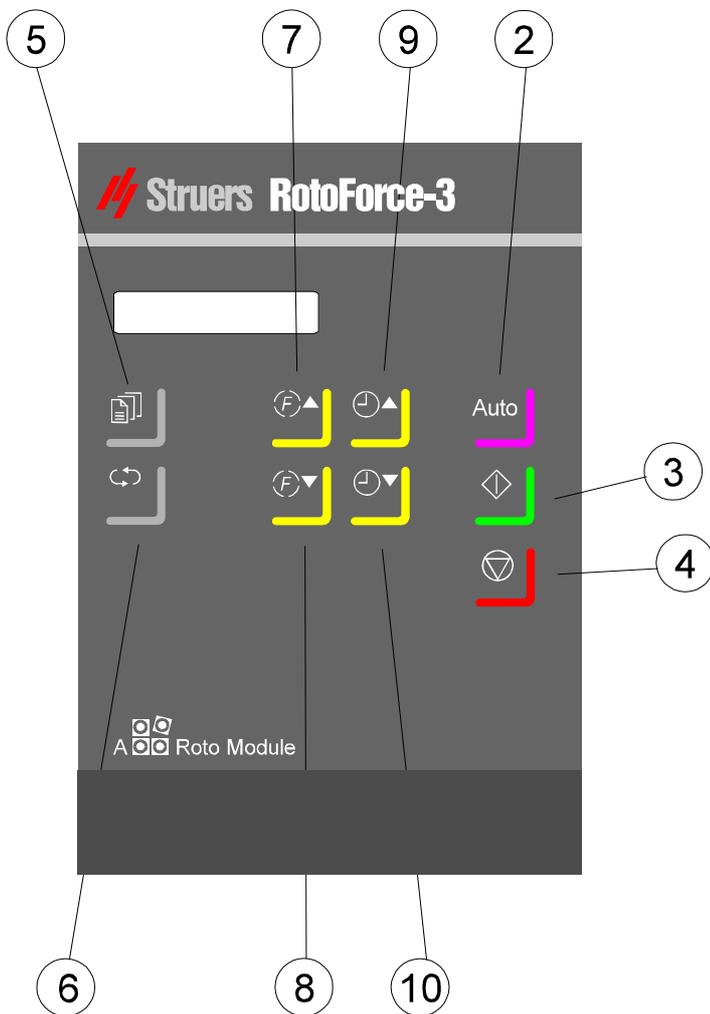
## **Mounting Lupo**

- Remove Lupo from its box
- Guide Lupo onto the connector on the supporting arm on RotoForce-3.
- Connect the nipple on Lupo to the nozzle underneath RotoForce-3 with the enclosed silicone tube.
- Remove the top lid of the bottle and fill with lubricant.
- Re-mount the top lid, and Lupo is ready for use.


## 2. Operation

Using the Controls  
Front Panel Controls of  
RotoForce-3 and Lupo

### RotoForce-3






### Lupo

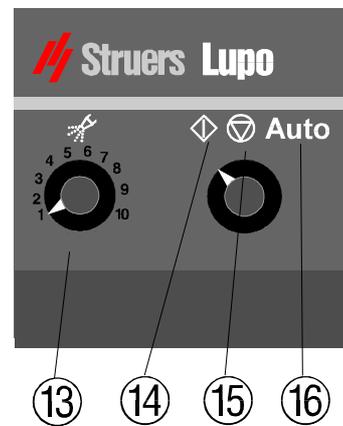
*RotoForce-3  
Instruction Manual*

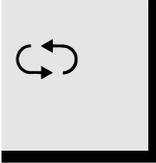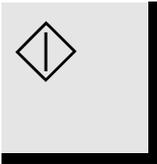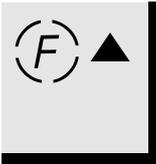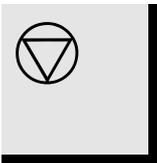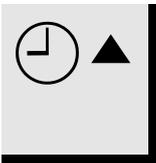**Front Panel Controls**  
*RotoForce-3*

Name	Key	Function	Name	Key	Function
① MAIN SWITCH		The main switch is located on the back of RotoForce-3. Turn clockwise to switch on the power.	⑤ MENU		Activates the set-up menu or moves down in the menu structure.
② AUTO		Turns Auto mode on or off (communication between the modules in the RotoSystem).	⑥ SELECT		Change the parameter of the installation menu.
③ START		Starts RotoForce-3.	⑦ ⑧ FORCE		Select FORCE: ▲ Increases the force ▼ Decreases the force
④ STOP		Stops RotoForce-3 Exits the set-up menu.	⑨ ⑩ TIME		Select TIME: ▲ Increases the time ▼ Decreases the time

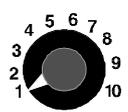*Lupo*

Name	Key	Function	Name	Key	Function
⑬ DOSING SWITCH		Switches between different step dosing levels.	⑭ START		Manual start of Lupo. When not in Auto mode Lupo can be used for manual pre-dosing.
(⑭, ⑮, ⑯) CONTROL SWITCH		Switches between Auto, Start and Stop mode. Lupo is supplied with power direct from RotoForce-3 and does not need a separate power switch.	⑮ STOP		Manual stop of Lupo. This mode does not stop neither RotoForce-3 nor the grinding/polishing machine.
			⑰ AUTO	<b>Auto</b>	When Auto mode is active start and stop of Lupo is controlled from RotoForce-3. When RotoForce-3 is started/stopped Lupo also starts/stops.

**Location of Main Switch**

① The main switch is located on the back of RotoForce-3.

**General Use of Keys**

- MENU , and SELECT  can only be activated when the machine has been stopped.
- SELECT  can only be activated when MENU  is selected.

*Acoustic Signals*

When pressing a key, a short beep indicates that the command has been accepted, whereas a long beep indicates that the key cannot be activated at the moment.

**The Handle**

The handle of RotoForce-3 is used to lock it in working position.

*Handle in Top Position:* when the handle is moved up into its top position, RotoForce-3 can be moved horizontally and placed correctly over the preparation disc.

*Handle in Bottom Position:* when RotoForce-3 is positioned correctly, move the handle down to lock RotoForce-3 in its working position.

**Changing to Another Lupo Unit (Accessory)**

- Set the mode switch on Lupo to STOP  position.
- Remove the current Lupo. Remove the silicon tube from the nipple underneath Lupo.
- Place the other Lupo (accessory) with an alternative lubricant in the connector on the supporting arm. Connect the silicone tube to the nipple.

**Changing Lubricant Type /  
Refilling**

- Set the mode switch on Lupo to STOP  position.
  - Remove Lupo from RotoForce-1.
  - Hold the Lupo lubricant bottle firmly and remove the top lid.
  - Empty the bottle, still mounted on Lupo, of any remaining lubricant.
    - Fill the bottle with mild, soapy water.
    - Put Lupo back on the supporting arm.
    - Set the mode switch on Lupo to START .
    - Set the dosing control to max. dosing (step 10) and let the water run through the nozzle for 2-3-min.
- Repeat the above points but fill with tap water to rinse out any soap residue.
- Refill with lubricant and remount the top lid.

***IMPORTANT***

Always avoid turning the socket cap while unscrewing the Lupo bottle, as the turning movement may cause damage to the socket cap gasket.

## Running a Preparation Method

### *Inserting a Specimen Holder*

- Place the specimen holder under the quick coupling.
- Press and hold the release button on the column of RotoForce-3, while guiding the pressure tap of the specimen holder into the coupling.
- Turn the specimen holder until the three pins engage in the corresponding holes.
- Release the button.

### *Adjusting the Position of RotoForce-3*

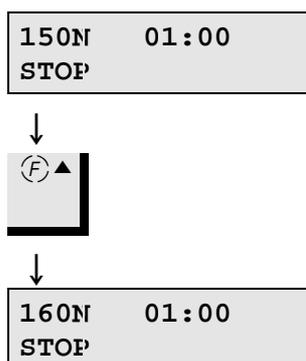
The position of the specimen holder must be adjusted correctly in relation to the preparation disc to achieve the best possible preparation results. This is done by moving RotoForce-3 horizontally.

- Generally the specimen holder should be positioned so that the specimens run close to the periphery of the preparation disc. A distance of about 2-4 mm is appropriate.
- There are three exceptions only: Grinding paper with double adhesive back, disc with magnetic backing and Petrodisc-M. In all three cases, the specimens should be protruding over the edge.


*NB:* In the case of Petrodisc-M, the specimens should protrude over the centre of the disc as well. This is done to achieve the best possible planeness of the specimen, and to maintain the planeness of Petrodisc-M

### *Adjusting the Force*

Press FORCE  $\text{\textcircled{F}}$ ▲ or  $\text{\textcircled{F}}$ ▼ to adjust the force to the desired value.

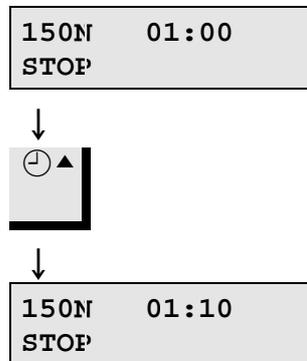*Adjusting the Preparation Time* Press TIME ⊖▲ or ⊖▼ to adjust the preparation time.

#### **Operation of Lupo**

- Adjust the dosing level
- Set Lupo to Auto mode

#### **Operation of RotoForce-3**

- After setting the correct time and force on RotoForce-3 and the dosing level on Lupo, press Auto on RotoForce-3.
- Press START ◊ on the RotoPol grinding/polishing machine to start both machines simultaneously.
- When the pre-set time has elapsed, all machines will stop simultaneously.

*RotoForce-3*  
*Instruction Manual*

# Reference Guide

Table of Contents	Page
<b>1. Force Reduction .....</b>	<b>16</b>
<b>2. Accessories.....</b>	<b>17</b>
<b>3. Trouble-shooting .....</b>	<b>19</b>
<b>4. Maintenance</b>	
Daily .....	20
Weekly.....	20
RotoForce .....	20
Lupo.....	20
<b>5. Technical Data.....</b>	<b>21</b>

## 1. Force Reduction

RotoForce-3 is equipped with the possibility of automatic reduction of the force towards the end of the preparation.

If FORCE REDUCTION is enabled, the force will automatically be reduced from the actual value down to 5 N during the last 30 seconds of the preparation step. This will result in lower material removal and smaller scratches than usual. The time for the subsequent preparation steps can be reduced as the amount of damaged material to be removed will be reduced.

1500N 04:00  
STOP

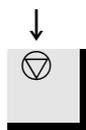
Press MENU  until you reach the FORCE REDUCTION screen.

FORCE REDUCTION  
NO

Press SELECT  to toggle between NO and YES.

FORCE REDUCTION  
YES

Press STOP  to accept the choice. Operation Menu will appear again.

150N 04:00  
STOP

## 2. Accessories

Specification	Code
<p><i>Specimen holder, ø 140 mm</i> for 6 specimens, ø 10-32 mm, aluminium for 6 specimens, ø 10-32 mm, stainless steel for 4 specimens, ø 10-40 mm, aluminium for 4 specimens, ø 10-40 mm, stainless steel</p>	<p>PEDAL PEDST PEDFI PEDIS</p>
<p><i>Specimen Holders ø 160 mm</i> for 6 specimens, each ø 25 mm/1", aluminium for 6 specimens, each ø 30 mm, aluminium for 6 specimens, each ø 1.1/4", aluminium for 4 specimens, each ø 40 mm, aluminium for 4 specimens, each ø 1.2", aluminium</p>	<p>MAXTO MAXUM MAXQU MAXLI MAXHA</p>
<p><i>Specimen Holders ø 160 mm</i> for 6 specimens, ø 12-40 mm, stainless steel for 6 specimens, ø 12-40 mm, aluminium for 12 specimens, ø 10-25.5 mm, stainless steel for 6 flat specimens (max. 34x60 mm), stainless steel for 6 specimens, ø30 or 22x44 mm, stainless steel for 3 flat specimens (max. 41x72 mm), stainless steel With 3 clamps for wire or plate (max. 24 x 52 mm) stainless steel</p>	<p>MAXCY MAXLU MAXDU MAXYK MAXIB MAXDI MAXAB</p>
<p><i>Specimen Holders ø 160 mm without Holes</i> plane, solid specimen holder for adhesion of specimens or for custom-made shapes Stainless steel Aluminium</p>	<p>MAXSO MAXAS</p>
<p><i>Specimen Holders ø 200</i> for 5 specimens, ø 27-55 mm, stainless steel for 3 specimens up to ø 70 mm or one specimen up to ø 180 mm, stainless steel for 3 flat specimens (max. 53 x 70 mm), stainless steel</p>	<p>MAXUK MAXRU MAXOT</p>
<p><i>Specimen Holders ø 200 mm without Holes</i> plane, solid specimen holder for adhesion of specimens or for custom-made shapes Stainless steel Aluminium</p>	<p>MAXON MAXAN</p>

*RotoForce-3*  
*Instruction Manual*

<b>Specification</b>	<b>Code</b>
<i>Uniforce, basic model</i> Device for levelling of specimens in specimen holders ø 140, 160, and 200 mm	MAXWI
<i>Uniforce, advanced model</i> With pressure foot for positioning the specimens	MAXFU
<i>Uniforce extension kit</i> Pressure foot for positioning the specimens	MAXMO
<i>Levelling Device for levelling of Specimens</i> In specimen holders 140 mm dia.	PEDFU
LUPO, optional extra drip lubricator	PEDNY

***Remember...***  
Struers offers a comprehensive range of consumables for both grinding and polishing. Please ask for separate leaflets.

### 3. Trouble-shooting

Display/Error	Cause	Action
<b>RotoForce-3</b>		
OVERLOAD	RotoForce-3 is stopped due to overloading of the motor.	Wait for the motor to cool down and the PAUSE/OVERLOAD message will appear on the display.
PAUSE/ OVERLOAD	The motor has cooled down after overloading and RotoForce-3 is ready for use.	Restart RotoForce-3.
PAUSE/ NO AIR	Air pressure too low.	Supply correct air pressure and restart.
LOCK HANDLE	The handle has not been pulled and RotoForce-3 is at the moment unlocked.	Lower handle.
Japanese Texts	RotoForce-3 has the possibility of showing the screen texts in Japanese. When choosing a European language in the LANGUAGE menu, you might accidentally press  MENU at the Japanese option.	To return to English, German or French texts, just follow this procedure: <ul style="list-style-type: none"> <li>■ Press STOP .</li> <li>■ Press MENU  two times.</li> <li>■ Press SELECT  until you reach the language you prefer.</li> <li>■ Press STOP .</li> </ul>
<b>Lupo</b>		
Bad or missing dosing	Dosing nozzle clogged.	Clean the nozzle with alcohol.
Lubricant leaking below the bottle	The bottle has not been fastened correctly	Fasten tightly. <b>Remember</b> to hold back the socket cap itself to avoid damage to the bottom gasket

## **4. Maintenance**

### **Daily**

Clean all accessible surfaces of the machine with a moist cloth

### **Weekly**

#### *RotoForce*

- Unlock RotoForce-3.
- Press START  twice.
- Clean the coupling with a clean cloth. An acid-free oil is recommended for the lubrication of the coupling to secure easy movement.
- Press STOP  to return the coupling into the cylinder.

#### *Lupo*

To avoid clogging of the Lupo nozzle, you should clean the nozzle once a week:

- Set the mode switch on Lupo to STOP  position.
- Remove Lupo from the supporting arm.
- Hold the Lupo lubricant bottle firmly and remove the top lid.
- Empty the bottle, still mounted on Lupo, of any remaining lubricant.
  - Fill the bottle with mild, soapy water.
  - Put Lupo back on the supporting arm.
  - Set the mode switch on Lupo to START .
  - Set the dosing control to max. dosing (step 10) and let the water run through the nozzle for 2-3-min.
  - Repeat the above points but fill with tap water to rinse out any soap residue.
  - Refill with lubricant and remount the top lid.

## 5. Technical Data

Subject			Specifications	
			Metric/International	US
Disc	Rotational speed		150 rpm	150 rpm
	Motor output 50/60Hz		90/80W	0.11/0.12HP
	Torque at disc	Continuous 50/60HZ	5.1/5.3Nm	3.8/3.9lb-ft
		Max. 50/60HZ	12.6/12.9Nm	9.3/9.5lb-ft
	Air pressure		6 bar	6 bar
Specimens	Force on specimen holder		30-400 N in 10N steps	30-400 N in 10N steps
	Lifting force, upwards		max. 60N	max. 60N
Dimensions and Weight	Width		270 mm	10.6"
	Depth		450 mm	17.7"
	Height		500 mm	19.7"
	Weight		25 kg	55.1 lbs

Subject		Specifications
Software and electronics	LC display	2x16 characters
	Controls	Touch pad
	Memory	EPROM, RAM and non-volatile RAM
	Automatic start/stop	With: Lupo
	Communication including automatic start/stop	With: RotoPol-21/22/25 Multidoser
	Data exchange	With: RotoCom
Environment	Safety Standards	General 89/392/EEC & 91/386/EEC & 93/44/EEC EN292-1, EN292-2, EN60204-1 (IEC 204-1) EMC 89/336/EEC & 92/231/EEC EN50081-1, EN50082-1 FCC part 15, subpart B
	Noise Level	Approx. 42 dB (A) at idle running, at a distance of 1.0 m/39.4" from the machine
	Surrounding temperature	5-40°C/41-104°F
	Humidity	0-95% RH non condensing

*RotoForce-3*  
*Instruction Manual*

<b>Subject</b>		<b>Specifications</b>
Supply Voltage	Power consumption 50/60HZ	180/165W (0.24/0.22HP)
	No. of phases	3 (3L + PE)
	Voltage / frequency:	Max. continuous load
	3 x 200 V, 50 Hz	1.35A
	3 x 220 V, 50 Hz	1.25A
	3 x 380 V, 50 Hz	0.80A
	3 x 415 V, 50 Hz	0.80A
	3 x 200 V, 60 Hz	1.40A
	3 x 220 V, 60 Hz	1.30A
3 x 440 V, 60 Hz	0.76A	
3 x 480 V, 60 Hz	0.72A	

# Quick Reference Card

## Inserting a Specimen Holder

- Place the specimen holder under the quick coupling.
- Press and hold the release button on the column of RotoForce-3, while guiding the top of the specimen holder into the coupling.
- Turn the specimen holder until the three pins engage in the corresponding holes.
- Release the button.

## Adjusting the Force

- Press FORCE  $\text{\textcircled{F}}^{\blacktriangle}$  or  $\text{\textcircled{F}}^{\blacktriangledown}$  to adjust the force to the desired value.

## Adjusting the Preparation Time

- Press TIME  $\text{\textcircled{T}}^{\blacktriangle}$  or  $\text{\textcircled{T}}^{\blacktriangledown}$  to adjust the time to the desired value.

## Adjusting the Position of RotoForce-3

- Unlock RotoForce-3 by moving the handle up.
- Adjust the RotoForce-3 so that the specimens are running at about 2 - 4 mm from the periphery of the preparation disc.
- Lock RotoForce-3 in that position by moving the handle down.

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Manual No.: 15257003

Date of Release 30.01.2001

*RotoForce-3/-4*  
*Spare Parts and Diagrams*

---

**Always state *Serial No* and *Voltage/frequency*  
if you have technical questions or when ordering spare parts.**

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**Instruction Manuals:** Struers Instruction Manuals may only be used in connection with Struers equipment covered by the Instruction Manual.

**Service Manuals:** Struers Service Manuals may only be used by a trained technician authorised by Struers. The Service Manual may only be used in connection with Struers equipment covered by the Service Manual.

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Some of the drawings may contain position numbers not used in connection with this manual.

*RotoForce-3/-4*  
*Spare Parts and Diagrams*

**Spare Parts List, RotoForce-3**

<b>Drawing</b>	<b>Pos.</b>	<b>Spare Part</b>	<b>Cat. no:</b>
14800107		<i>Rotating head, complete</i>	
	20	O-ring 68.00-2.00 (10 pcs.)	RIO20080
	30	Key + half moon ring	R4790418
	40	Pressure spring Ø22.1x0.9	R4790421
	110	K-ring N 896/55.1	RIK00055
	130	Rubber pad (10 pcs)	R4790442
15240061		<i>Front Cabinet, complete</i>	
	30	PCB, with PROM & Display	15242398
	30	PROM, latest version	R5253901
	30	Clamping ring for transducer on PCB	14590110
15240051		<i>Back Cover, complete</i>	
	110	Main switch, complete	2SE20401
	380	630mA Fuse (10 pcs)	315MP138
15240031		<i>Chassis, complete</i>	
	80	Ball bearing 6007-2z ø35/62x14	2BK00112
	140	Set of seals	14790103
	240	Cover Ø 21 (10 pcs)	RGD54025
	310	Poly-V belt (L=508 mm)	RJD30508
	400	Solenoid valve block (Y4-Y5), complete	15240032
	400	3/2 Solenoid valve 12VDC	RYM10114
	400	Gasket, PVC M5 (10 pcs)	14792902
430	Micro switch	14790217	

*RotoForce-3/-4*  
*Spare Parts and Diagrams*

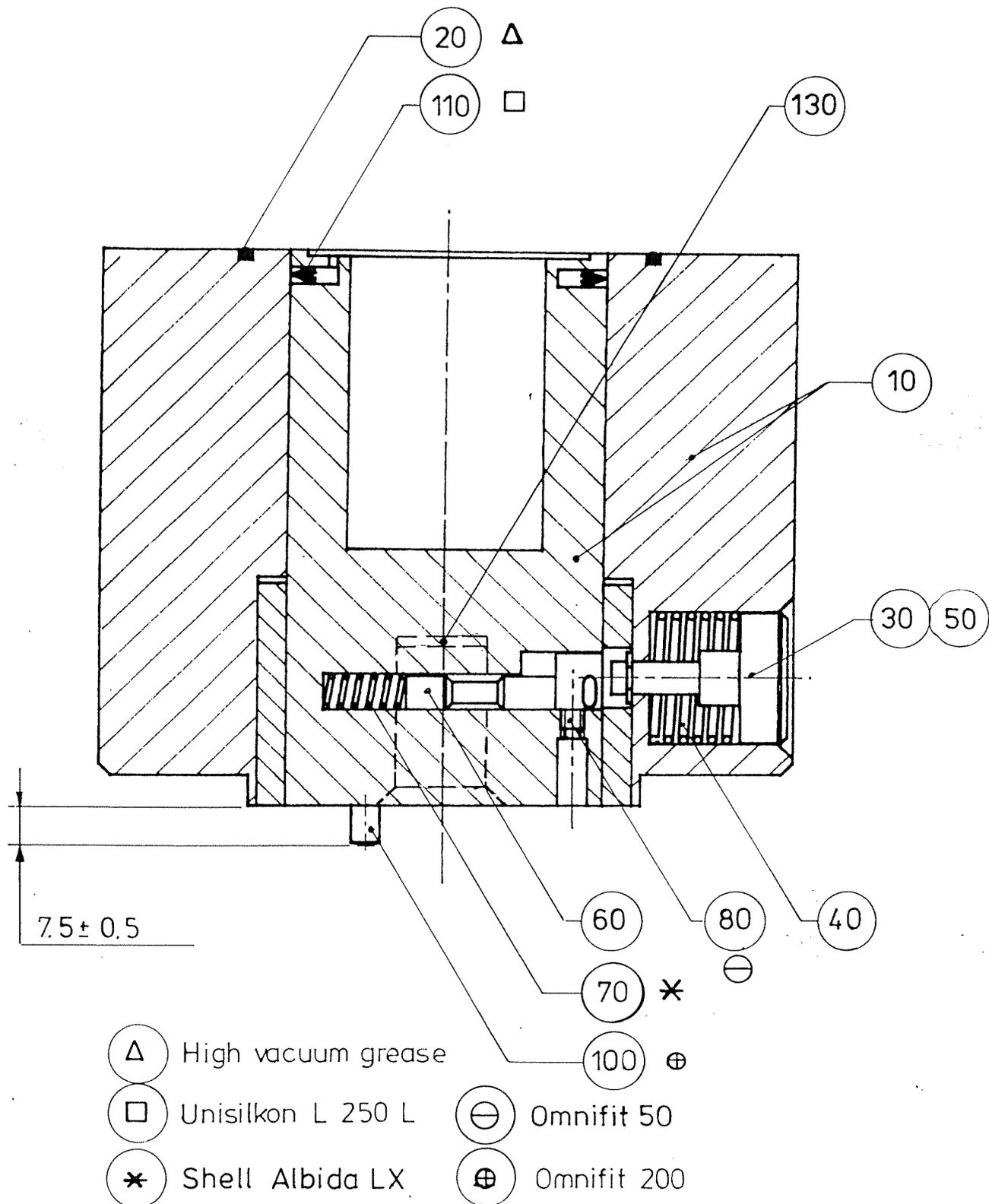
**Spare Parts List, RotoForce-4**

<b>Drawing</b>	<b>Pos.</b>	<b>Spare Part</b>	<b>Cat. no:</b>
15250032		<i>Chassis, complete</i>	
	60	Ball bearing, 6007-2Z	2BK00112
	120	Set of seals	14790103
	a	O-Ring 15.60-2.40 (2 pcs)	
	b	O-Ring 7.65-1.78 (2 pcs)	
	c	O-Ring 28.00-2.00	
	260	Cover Ø 21 ELMA (10 pcs)	RGD54025
	370	Poly-V belt (L=508 mm)	RJD30508
	420	V-ring VS-0040	RIV00040
	430	Set of seals	14790103
	450	QUAD-Ring	RIQ04215
	545	Gasket, PVC M5. (10 pcs.)	14792902
	550	3/2 Solenoid valve M5 12VDC	RYM10112
	570	3/2 Solenoid valve 12VDC	RYM10114
	610	Micro switch	14790217
	740-750	Screw with ball and spring, M8 + Nut	RTX90816
	14790390		Rotating head, complete
		Replacement Rotating head, complete, RF-4	14792904
20		O-Ring 110.00-2.00	RIO20090
30		O-Ring 36.00-2.00 (5 pcs)	386MP025
40		Key + half moon ring	R4790418
140		Rubber pad (10 pcs)	R4790442
180		Bellows, rubber/V6-752	454MP078
		Clamping ring for foot (6 pcs)	R3750351
210		Pressure foot	454MP077
220		O-ring Ø8.30-2.40	RI024015
15250062			<i>Front Cabinet, complete</i>
	10	Front cabinet with touch pad keyboard	15250500
	30	PCB with PROM & Display	15252372
	30	PROM, latest version	R5253901
15250052		<i>Back Cover, complete</i>	
	70	Mechanical blocking CM 4	2KH10021
	80	Main switch, complete	2SE20401
	250	630mA Fuse (10 pcs)	315MP138
14790020		<i>Pneumatic connections</i>	
	10	Quick coupling double	2NF10055
	11	Spiral hose type SPF 15, 333 mm	2WK90015


*RotoForce-3/-4*  
*Spare Parts and Diagrams*

**Common Parts to RotoForce-3 and RotoForce-4**


<b>Drawing</b>	<b>Pos.</b>	<b>Spare Part</b>	<b>Cat. no:</b>
14600230		<i>Fixture, complete</i>	
	20	Glider	14600232
	60	O-ring 3.00-1.00 (10 pcs.)	315MP036
	70	Nozzle fixture	R4600083
14820010		<i>Lupo</i>	
	30	Valve block	R4820021
	40	Nylon washer (20 pcs.)	R4822901
	50	Bottle screw	R4820026
	60	Washer Ø31x1 (10 pcs)	R4820027
	70	Lupo lubricant bottle	R4820028
	71	Top Lid for Lupo Bottle	R4820056
	80	Solenoid valve, 24VDC	RYM10124
	80	PCB logical card with solenoid valve	R4823000
	100-120	ELMA dial (mode switch) button, ø3	255MP074A
130	Dosing switch button	2GD12225	


Matr.		Målestok: 1:1					
		Ikke ang. tolerancer: ±	Overfl.-behandl. Ja <input type="checkbox"/> Nej <input checked="" type="checkbox"/>				
		Ikke ang. ruheder:	<table border="1"> <tr> <td>Date</td> <td>Sign.</td> </tr> <tr> <td>Tegn. 31.7.92</td> <td>B</td> </tr> <tr> <td>Kontr.</td> <td>O: B</td> </tr> </table>	Date	Sign.	Tegn. 31.7.92	B
Date	Sign.						
Tegn. 31.7.92	B						
Kontr.	O: B						
B: 7 ± 0,5 → 7,5 ± 0,5 Fedt → Shell Albida LX 00997 BMS/791	Kode	Anv. i type:	Erst.:				
A Omnifit 50 tilf. 16.11.92 B3/791		Hoved, saml.	14800107 E				

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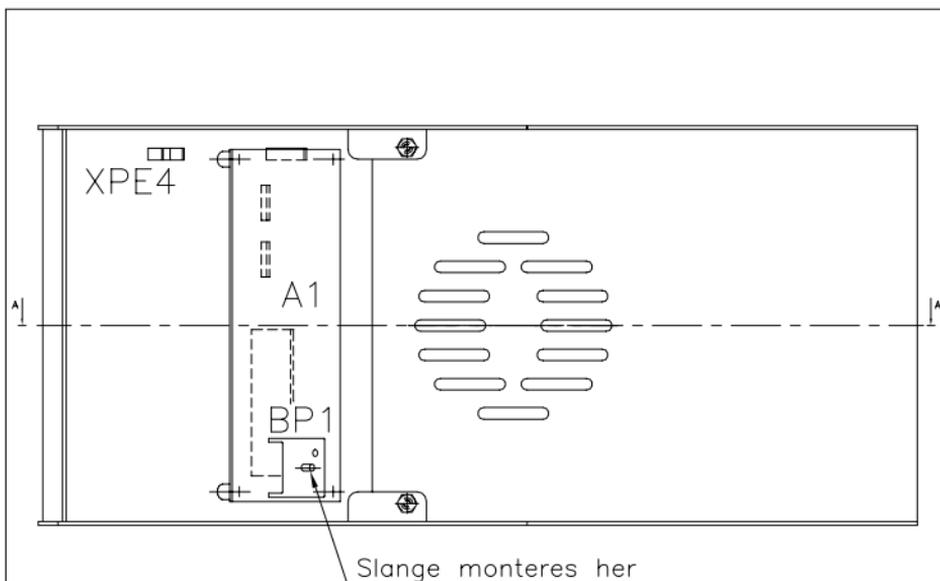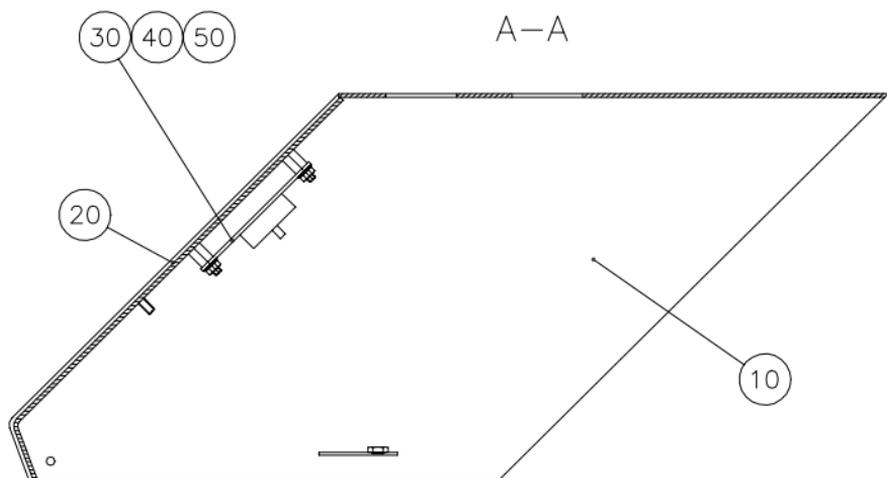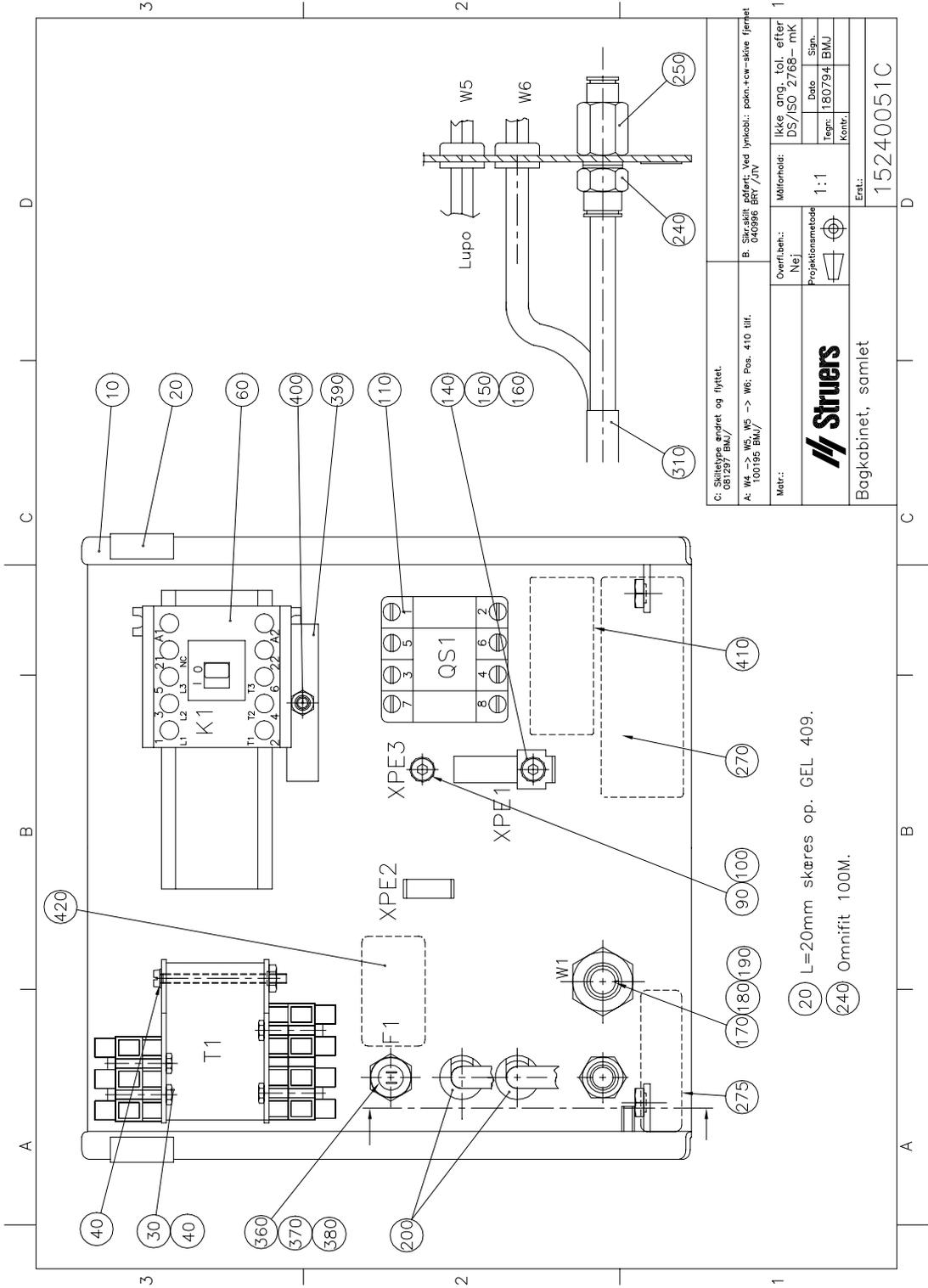
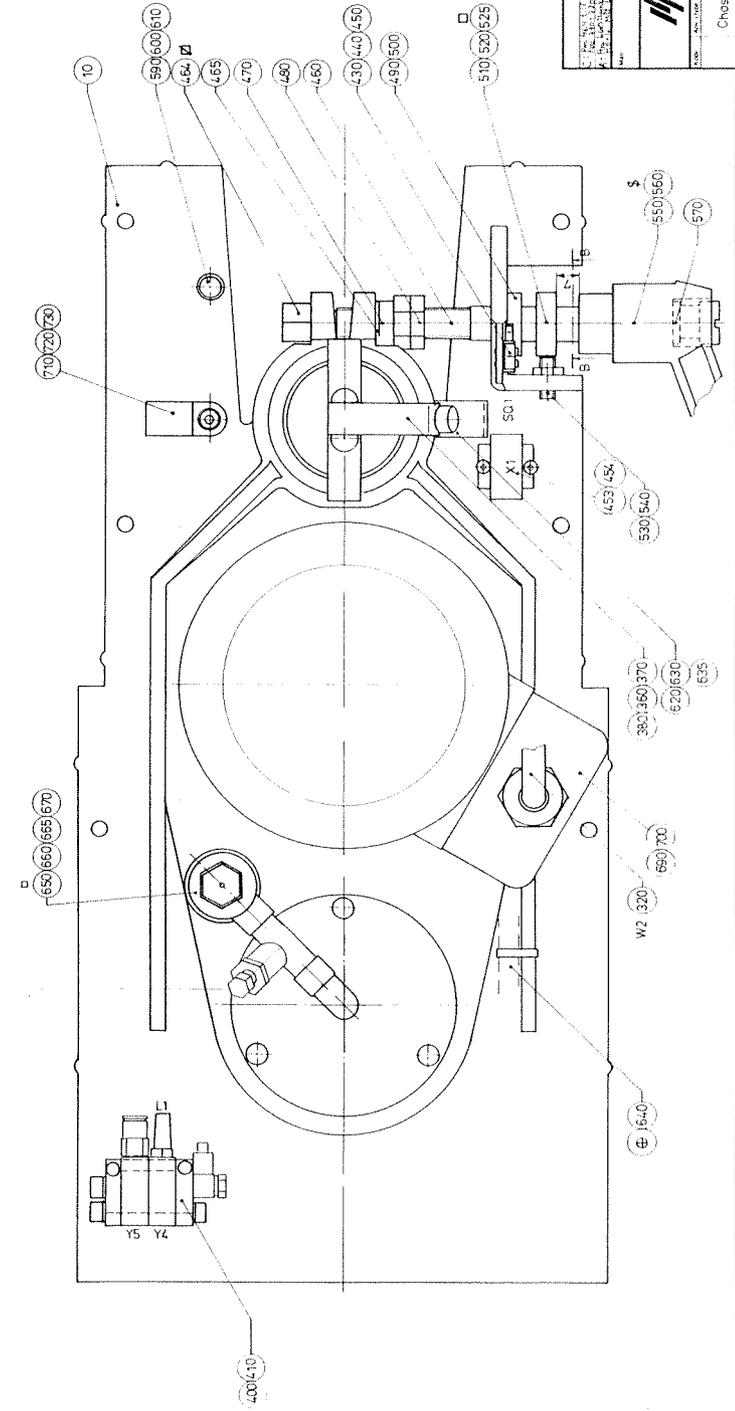
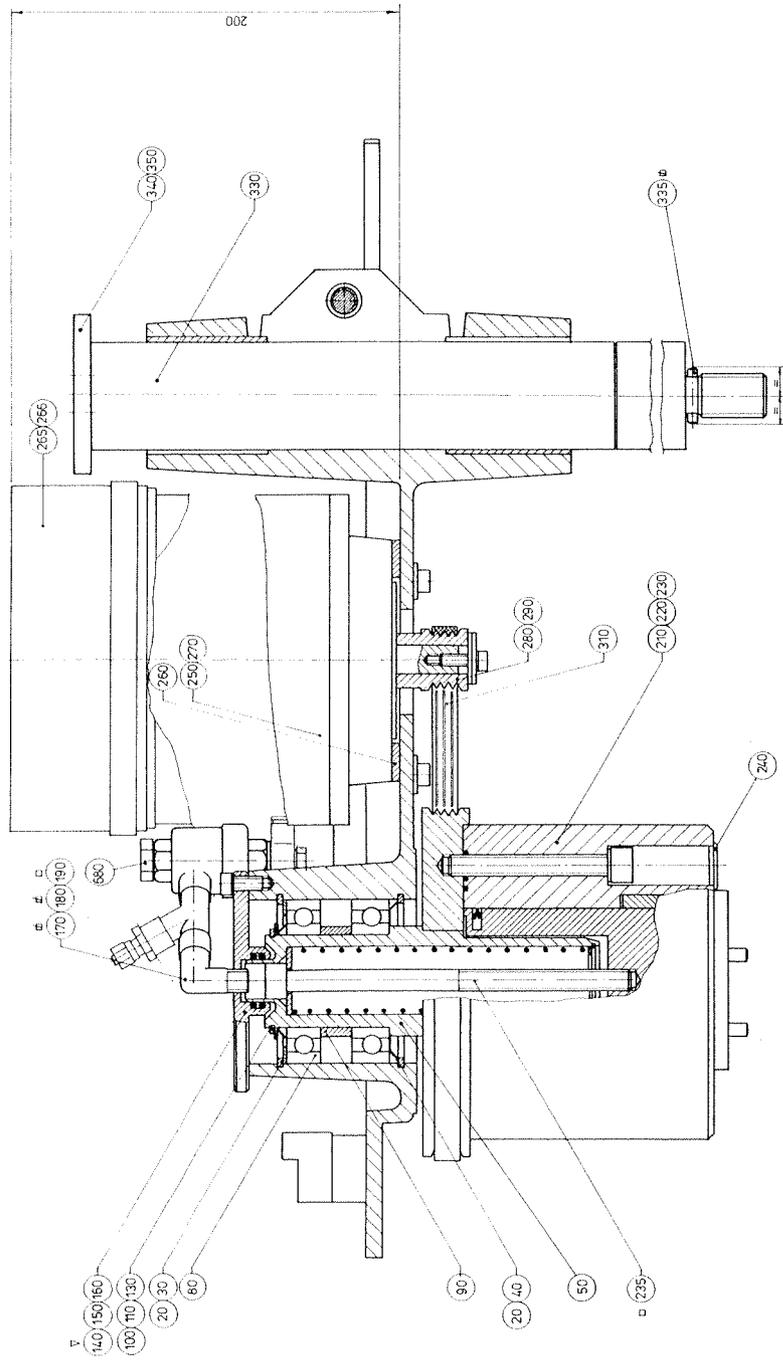
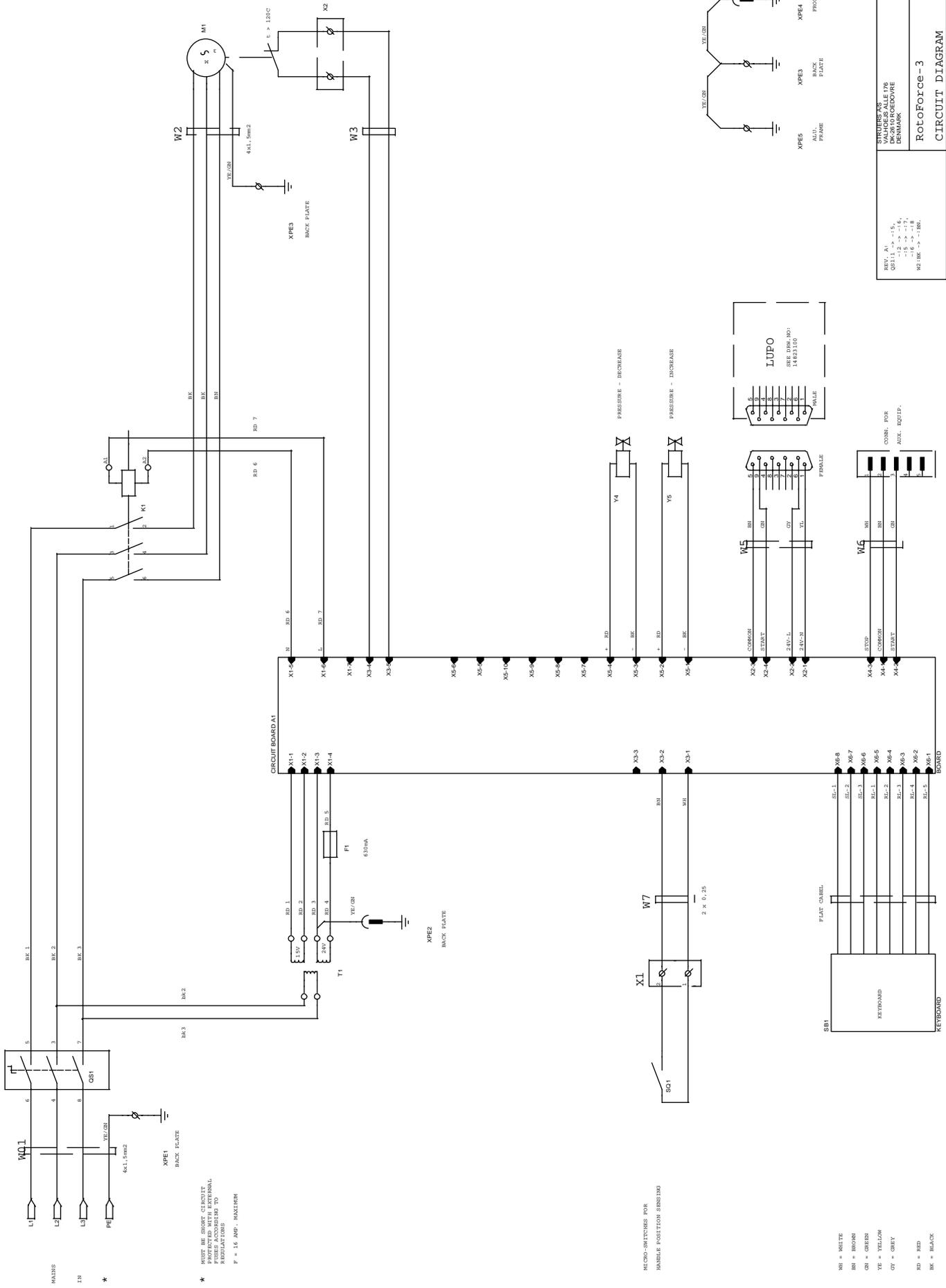
\* MUST BE SHORT CIRCUIT PROTECTED WITH EXTERNAL REGULATORS  
 P = 16 AMP. MAXIMUM

MICRO-SWITCHES FOR HANDLE POSITION SENSING  
 WH = WHITE  
 BR = BROWN  
 GN = GREEN  
 YE = YELLOW  
 OY = GREY  
 BK = RED  
 BK = BLACK

SYNERGY'S VALBORES ALLE 796 28000 COPENHAGEN DENMARK		<b>RotoForce-3</b> CIRCUIT DIAGRAM	
REV. A1 - 15	Scale	15243100	Rev
REV. A2 - 16	C	1	of 1
REV. A3 - 17	Sheet	1	
REV. A4 - 18	DWG NO	15243100	
REV. A5 - 19	Thursday, June 08, 2000		
REV. A6 - 20			

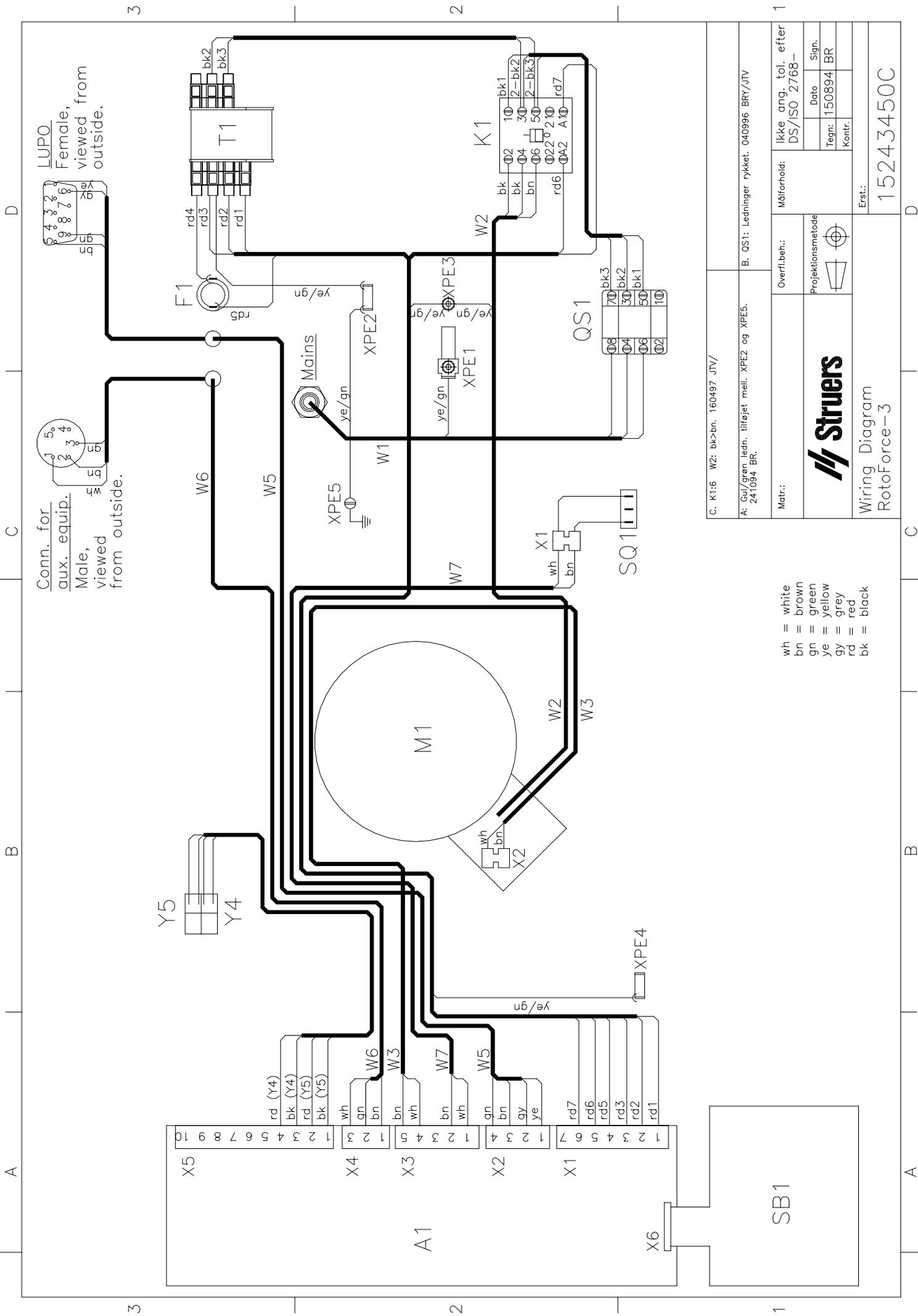Conn. for  
aux. equip.  
Male,  
viewed  
from outside.

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Herewith declares that

<i>Product Name:</i>	RotoForce-3
<i>Type No:</i>	524
<i>Machine Type:</i>	Specimen mover

is in conformity with the provisions of the following directives:

**Safety of Machinery** 89/392/EEC and 91/368/EEC and 93/44/EEC according to the following standard(s):  
EN292-1:1991, EN292-2:1991, EN60204-1:1992.

**EMC-Directive** 89/336/EEC and 92/31/EEC according to the following standard(s):  
EN50081-1:1991, EN50082-1:1992.

**Low Voltage Directive** 73/23/EEC and 93/68/EEC according to the following standard(s):  
EN60204-1:1992.

**Supplementary Information** The equipment complies with the American standards:  
FCC part 15, subpart B and UL508.

The above has been declared according to the global method, module A

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Klaus Kisbøll, Entwicklungsleiter, Struers A/S
Klaus Kisbøll, Chef du service developpement, Struers A/S